


High-Throughput Robot for Crystallization in Lipid Mesophase

Manufacturer: SIAS, AG, modified and assembled in the laboratory of Professor Martin Caffrey (see Ref)

Primary Use: Carrying out protein crystallization experiments using lipidic mesophases.

Description:

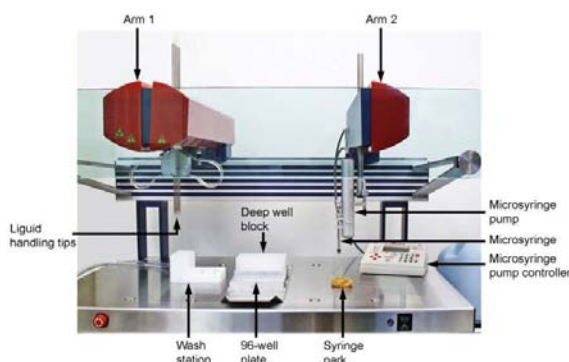

The robot was designed and built first and foremost to perform automated crystallization of membrane proteins using the *in meso* method. However, it also works with other crystallization protocols. The robot was built based on Xantus, a commercially available liquid handling robot (Sias AG, Switzerland; Zinsser NA, Northridge, CA, USA). The open style, modular design and software flexibility of the Xantus allowed its adaptation to perform the steps needed for *in meso* crystallization. The standard Xantus robot consists of a 1 X 0.7 m deck with two independent arms. Arm 1 is used for liquid aspiration and dispensing, while arm 2 is typically a plate gripper and moving

device. Arm 1 is equipped with four small-volume liquid-handling tips for pick-up and placement of precipitant solutions. Arm 2 handles the viscous protein/lipid mesophase. To this end, the gripper device uses a motor-driven positive-displacement Hamilton syringe (UMP II with Micro4 controller, World Precision Instruments, Sarasota, FL, USA). The controller for the latter was interfaced to the Xantus through an RS-232 cable. The positional accuracy of arm 2 was raised by increasing the gear ratio of the z motor by a factor of 8.

This LCP robot has been installed in the JCIMPT laboratory of Professor Ray Stevens on long-term loan from Ohio State University as part of collaboration with Professor Caffrey's laboratory. This robot is used routinely for preparing crystallization experiments on GPCR protein detergent complexes; it was used in producing high quality crystals of the GPCR's β 2-adrenergic receptor and Adenosine 2A receptor resulting in the determination of their 3-dimensional structure.

This robot is currently not available commercially. Those interested in building a system can contact Sias, AG. JCIMPT is working with Sias, AG and Zinsser NA to design and manufacture a new production version of the robot.

References

1. Cherezov, V., Peddi, A., Muthusubramaniam, L., Zheng, Y.F. & Caffrey, M. A robotic system for crystallizing membrane and soluble proteins in lipidic mesophases. *Acta Crystallogr D Biol Crystallogr* **60**, 1795-807 (2004)
2. Cherezov, V., and Caffrey, M. (2007). Miniaturization and automation for high-throughput membrane protein crystallization in lipidic mesophases. In: Protein crystallization strategies for structural genomics. N.E. Chayen, editor. International University Line: San Diego.



JCIMPT

JOINT CENTER FOR INNOVATIVE MEMBRANE PROTEIN TECHNOLOGIES

3. Cherezov V., Rosenbaum D.M., Hanson M.A., Rasmussen S.G., Thian FS, Kobilka TS, Choi HJ, Kuhn P, Weis WI, Kobilka BK, Stevens RC., High-resolution crystal structure of an engineered human beta2-adrenergic G protein-coupled receptor. *Science*. (2007) **318**(5854):1258-65
4. Hanson MA, Cherezov V, Griffith MT, Roth CB, Jaakola VP, Chien EY, Velasquez J, Kuhn P, Stevens RC.. "A specific cholesterol binding site is established by the 2.8 Å structure of the human beta2-adrenergic receptor.", *Structure*. 2008Jun;**16**:897-905.
5. Jaakola VP, Griffith MT, Hanson MA, Cherezov V, Chien EY, Lane JR, Ijzerman AP, Stevens RC., "The 2.6 angstrom crystal structure of a human A2A adenosine receptor bound to an antagonist.", *Science*. (2008);**322**:1211-7